

X-ray crystal spectrometer

M. Goto, T. Oishi, Y. Kawamoto

e-mail: goto@nifs.ac.jp

1. Objective

- [1] Ion temperature, $T_i(0)$
- [2] Toroidal rotation, $V_t(0)$

2. Apparatuses

2.1. Crystal holder

- Four curved crystals are installed on the crystal holder (see Fig. 1).

Quartz (2020): He-like ArXVII (3.95Å)

Quartz (2023): He-like TiXXI (2.61Å)

Quartz (3140): He-like CrXXIII (2.18Å)

Quartz (2243): He-like FeXXV (1.85Å)

- $T_i(0)$ is generally measured from ArXVII with Ar gas puffing

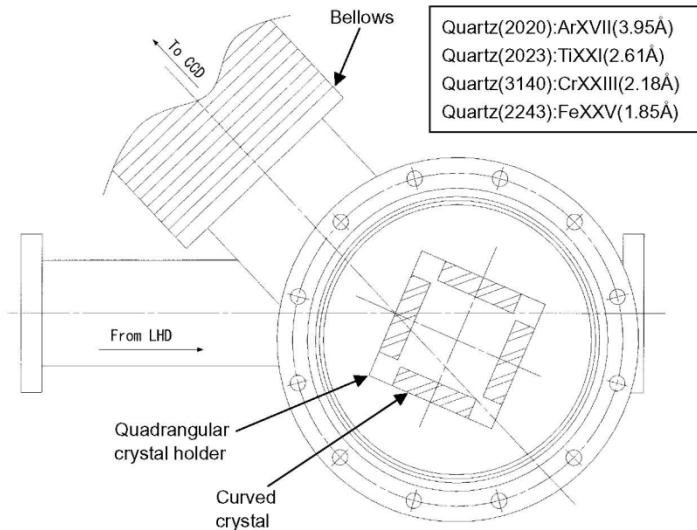


Fig. 1. Viewing chords of toroidal array

2.2. Port assembly

- Crystal spectrometer is installed on #3-O (Fig. 2)
- Observation area: equator plane ($-10\text{cm} \leq \Delta Z \leq +10\text{cm}$, ΔR depends on T_e profile)

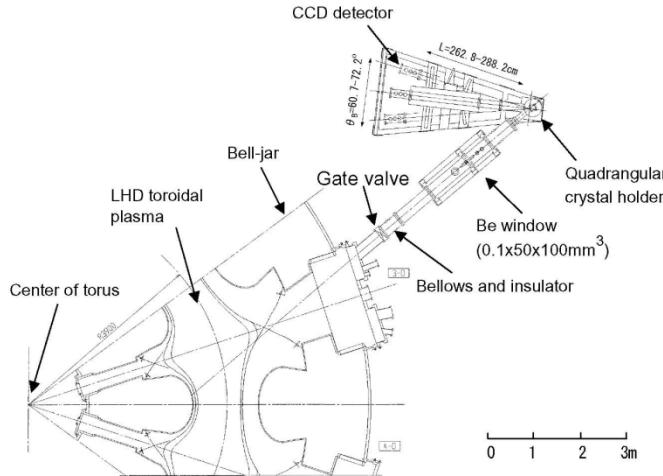


Fig. 2. Viewing chords at #5-O port

2.3. Specifications

- Time resolution: $\Delta t=5\text{ms}$

3. Operation of crystal spectrometer

- Before experiment: opening gate valve, turning on power supply and setting detector parameters
- After experiment: closing gate valve and turning off power supply
- Ar puffing is necessary at beginning of discharges

4. Available data by “Retrieve”

Kaiseki-data server

-ti_crystal

5. Remarks

- Example of Ti(0) measurement (see Fig. 3)

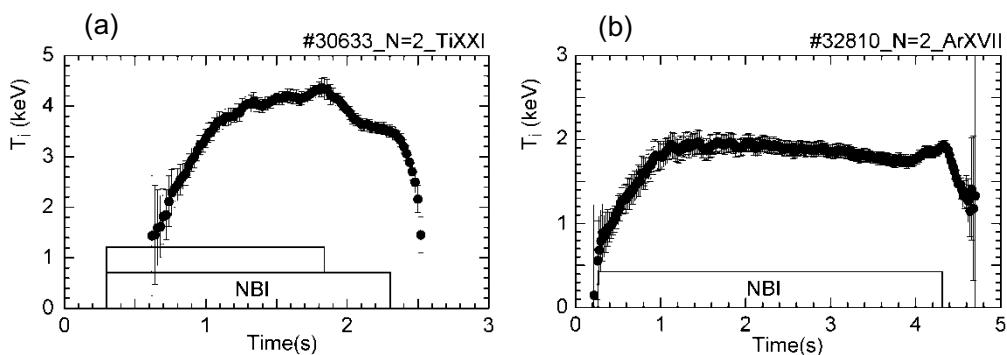


Fig. 3. $T_i(0)$ during NBI discharges for (a) TiXXI and (b) ArXVII

- Example of Doppler profiles of ArXVII during high-Ti discharges (see Fig. 4)

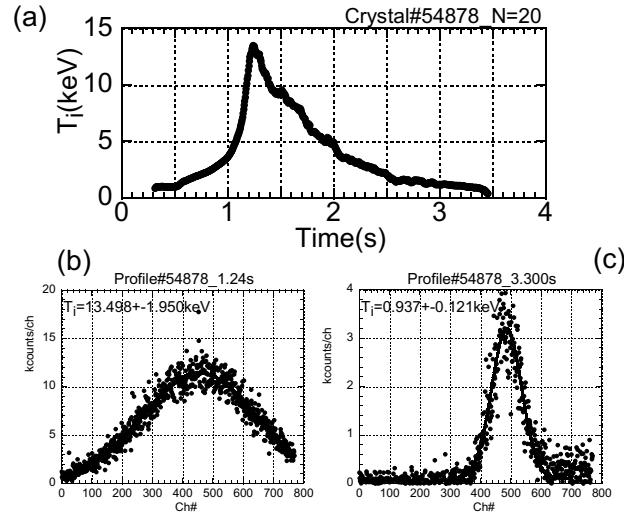


Fig. 4. Doppler-broadened ArXVII profiles; (a) $T_i(t)$ and ArXVII line profiles at (a) $t=1.24s$ and (c) $t=3.3s$.

- Example of He- and Li-like Ar spectrum (see Fig. 5)

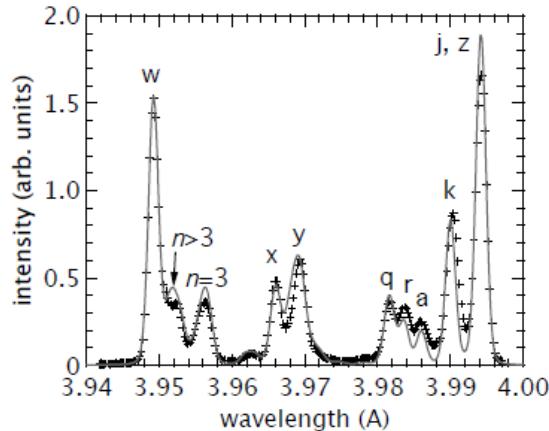


Fig.5. Typical example of He- and Li-like Ar spectrum.

References

- S.Morita, M.Goto et al., NF **42** (2002) 1179.
- S.Morita and M.Goto RSI **74** (2003) 2375.
- S.Morita, M.Goto et al., NF **43** (2003) 899.
- I.Sakurai, Y.Tawara et al., RSI **77** (2006) 10F328.
- I.Sakurai, Y.Tawara et al., PFR **2** (2007) S1068.
- M.Goto and S.Morita PFR **5** (2010) S1040.
- and others